


Caution: To ensure correct and safe use of the products, make sure that you also read the Operating Instructions on DVD carefully prior to operation.



How to clean the **OactiCAP**


1

After disconnecting all the electrodes from the cap, soak the cap in warm water mixed with a dime sized amount of baby shampoo.

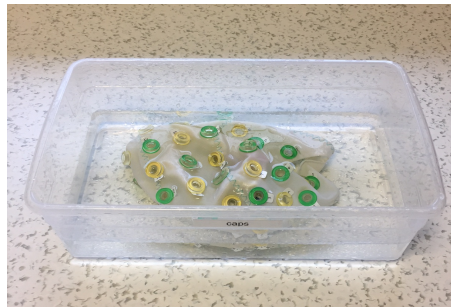

2

Wash the cap carefully, making sure that no gel remains on the cap

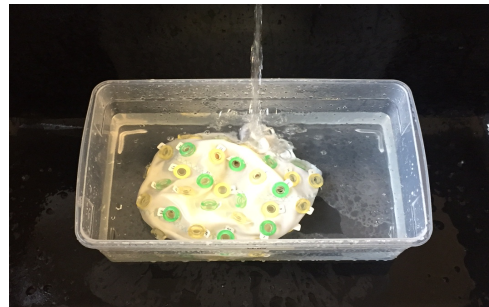

3

Rinse the cap carefully. Make sure no soap remains on the cap. The plastic rings may fall off. Make sure they don't go down the drain.



4

Put the cap on a paper towel, and lightly pat with another paper towel to dry.



5

Leave the cap on manikin to dry. If the cap is not dry when you want to use it, blow dry with low setting before using it.



Report to Lab Manager if:

- You lost a plastic ring on the cap.
- The cap is no longer fitting well.
- If the cap comes into contact with any potential infectious agent (for example, blood, contact the lab manager for disinfection procedure).